



Sterile Sample Bag (For Liquid)

【Application】

Used for aseptic collection of liquid samples (milk, soy milk, juice, purified water, mineral water, etc.).

【Advantage】

Ready to use after sterilization by gamma irradiation. No need to autoclave, wash and dry.

【Instructions】

Take a sterile water sample collection bag, unscrew the lid, and add the collected Liquid sample to the 200ml or 500ml **graduation line** of the bag (for the sample volume standard.) Tighten the lid.

【Sterilization Method】

Gamma-ray radiation sterilization

【Specifications】

AL-BG500E: 500ml, 10pcs/bag, 10bags/carton.

AL-BG200E: 200ml, 10pcs/bag, 10bags/carton.

【Shelf life】

Two years

【Precautions】

1. Hand pinch the neck of the bag mouth.
2. Gently shake the bag to be fully opened.
3. The line of sight is parallel to the 200ml or 500ml graduation scale.
4. Avoid contaminating the bag mouth with fingers or other objects.